



All-in-One™ miRNA qPCR Primer Validation Report

GeneCopoeia, Inc.
9620 Medical Center Drive, #101
Rockville, MD 20850
USA

301-762-0888
866-360-9531

inquiry@genecopoeia.com

www.genecopoeia.com

All-in-One™ miRNA qPCR Primers Sample Validation Report

- I. Materials and Methods
- II. Procedure
- III. Test Results
- IV. Limited Warranty

I. Materials and Methods

1. Instrument
iQ5 Real Time PCR Detection System: Bio-Rad

2. Reagents

All-in-One™ miRNA qRT-PCR Detection Kit (Catalog Nos. AOMD-Q020 or AOMD-Q050)
 Validation templates: cDNAs extracted from 10 different human tissues
 Detection primers: Total 6 pairs

Catalog number	Primer ID	Mature_Acc	Mature_ID	PCR size
miRQP0227	hsmq-0031	MIMAT0000069	hsa-miR-16	75bp
miRQP0074	hsmq-0032	MIMAT0000422	hsa-miR-124	73bp
miRQP0096	hsmq-0034	MIMAT0000423	hsa-miR-125b	75bp
miRQP0175	hsmq-0035	MIMAT0000429	hsa-miR-137	76bp
miRQP9901	hsnRNA-U6	NR_002752.1	RNU6B	80bp
MiRQP0293	hsmq-0655	MIMAT0000073	hsa-miR-19a	72bp

II. Procedure

RNA Extraction

1. Sample preparation

Place a small amount of material from a human tissue sample into a pre-cooled mortar. Add a small amount of liquid nitrogen and grind the tissue to a fine powder. Transfer the powder to a centrifuge tube containing 1 ml of TRIzol (Invitrogen). Vortex for 5 minutes. (If the samples are cultured cells, use about $10^6 \sim 10^7$ cells with 1 ml of TRIzol). Pipette up and down until the cells are completely lysed).

2. Phase separation

Leave the cell or tissue samples at room temperature for about 10 minutes. Add 200 μ l of chloroform per 1 ml of TRIzol. Close the cap and vortex vigorously for 1 minute. Let the samples settle down at room temperature for 2-5 minutes. Centrifuge at 12,000 g for 15 minutes at 6°C. Remove tubes from the centrifuge being careful not to disturb the liquid. The samples should be separated into 3 layers with RNA in the top layer.

3. RNA Precipitation

Carefully transfer about 450 μ l (per 1 ml TRIzol) of the supernatant to a new centrifuge tube containing 600 μ l of cold 2-propanol. Mix well and keep at -20°C for 10 minutes. Then centrifuge at 12,000 g for 10 minutes (at 6°C).

All-in-One™ miRNA qPCR Primer Validation Report

4. Washing

Remove the supernatant. Add 500 μ l of cold 75% ethanol. Vortex until a pellet forms at the bottom of the tube. Centrifuge at 12,000 g for 5 minutes (6°C). Remove the supernatant. Centrifuge briefly again and remove the remaining supernatant.

5. Dissolve RNA

Air-dry the samples for 5 to 10 minutes. (Do not over dry the samples. The samples are dry when they turn white). Dissolve the samples in 30 μ l of DEPC water. Label the samples properly and store them at -80°C.

6. Determine RNA Concentration

Take 1 μ l of RNA sample. Dilute the sample 1:10 in DEPC water. RNA concentration can be measured with a NanoDrop (Thermo Scientific). DEPC water should be used as a blank. Record both the RNA concentration and the ratio of A_{260}/A_{280} .

7. RNA electrophoresis

7.1. Prepare denaturing gel

Add 1 g of agarose into 75 ml of de-ionized water. Boil the agarose for 1–2 minutes and cool it down to about 70°C. Add 10 ml of 10X MOPS, 15 ml of formaldehyde and 5 μ l of ethidium bromide (EB). Pour the gel into a tray with big combs. Cover the tray.

7.2. Prepare electrophoresis buffer (1X MOPS)

Take 50 ml of 10X MOPS and dilute 1:10 to the final volume of 500 ml with de-ionized water. Pour the buffer into a gel box, and add some EB to it.

7.3. Preparation of RNA sample


Take 3 μ l of the RNA sample and add DEPC to a total volume of 18 μ l. Heat the sample at 65°C for 10 minutes. Cool it down immediately. Add 2 μ l of 1X RNA loading buffer into the sample.

7.4. RNA electrophoresis

Place the RNA gel in a gel box with 1X MOPS electrophoresis buffer. Pre-run the gel at 100 V for 5 minutes. Load the treated RNA samples into the gel and run at 100 V until the bromophenol blue runs to about one third the length of the gel. Take a picture using a UV scanner.

8. RNA detection results

Electrophoresis results of RNAs from 10 different tissues (3 μ l RNA each well). See table 8.2 for the specific lane information.

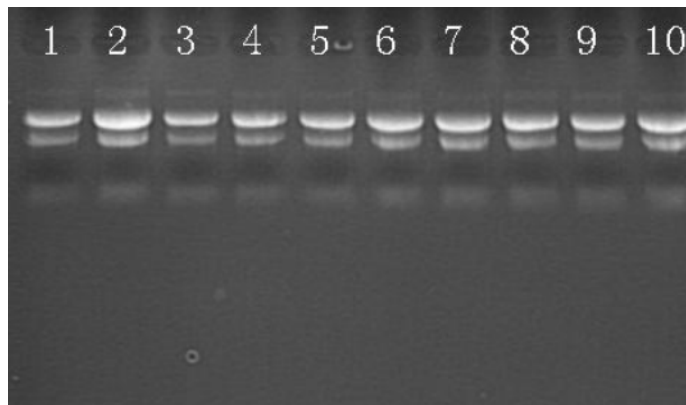

All-in-One™ miRNA qPCR Primer Validation Report

Table 8.2. RNA sample source, concentration and the ratio of OD260/OD280.

Lane	Tissue	Concentration (ng/μl)	OD A ₂₆₀ /A ₂₈₀
1	Brain	2385	1.9
2	Lung	2430	1.94
3	Liver	2521	1.91
4	Kidney	3444	1.94
5	Breast	2785	1.87
6	Testis	2972	1.9
7	Placenta	3515	1.91
8	Spleen	3344	1.91
9	Heart	3394	1.91
10	Pancreas	3101	1.91

Note: In order to detect miRNAs, the extracted RNA must contain small molecular weight RNAs. The kit used for RNA extraction must be applicable for isolation of total RNA or small RNA. The quality of RNA is critical to the success of downstream experiments. Follow the instructions of the RNA extraction kit exactly. Examine RNA samples by electrophoresis to ensure high quality.

miRNA reverse transcription

1. Thaw all the reagents needed for miRNA reverse transcription. Mix reagents well by gently inverting the tubes. Spin down briefly and keep on ice.
2. Prepare miRNA reverse transcription reaction: Add the following reagents into an RNase-free reaction tube which is pre-cooled on ice. The final volume should be 25 μl.

Reagents	Volume	Final concentration
Total RNA		2 μg
2.5 U/μl Poly A Polymerase	1 μl	2.5 U
RTase Mix	1 μl	
5X Reaction Buffer	5 μl	1X
ddH ₂ O (RNase/DNase free)	to 25 μl	

Note: The total RNA in the reaction must contain small molecular RNA. If total RNA is used, the amount of RNA should be between 1 ng ~ 5 μg. If the purified small molecular RNA is used, the amount of RNA should be between 0.1 ng ~ 1 μg

All-in-One™ miRNA qPCR Primer Validation Report

- Reverse Transcription Reaction: Mix reaction solution well. Spin down briefly. Incubate the reaction solution at 37°C for 60 minutes. Terminate the reaction by heating at 85°C for 5 minutes. The products of reverse transcription can be diluted 1:5 with sterile water for the downstream qPCR reaction.

qPCR to detect miRNA

- Thaw 2X All-in-One qPCR mix from the All-in-One miRNA qRT-PCR Detection Kit. Mix well by inverting the tube several times. Spin down briefly and keep on ice.
- Prepare qPCR reaction solution on ice. All miRNAs are tested in duplicate. NTC (No template control) is tested singly.

Reagents	Volume	Final Concentration
2×All-in-One qPCR Mix	10 µL	1X
All-in-One miRNA qPCR Primer	(2 µM) 2 µL	0.2 µM
Universal Adaptor PCR Primer	(2 µM) 2 µL	0.2 µM
First strand cDNA (diluted 1:5)	2 µL	
ddH ₂ O	4 µL	
Final volume	20 µL	

- Mix the qPCR reaction solution well. Transfer the solution to a PCR tube. Spin briefly to make sure that the reaction solution is at the bottom of the tube.
- The qPCR reaction, using the standard 3-step method to detect the DNA amount (these experiments were designed based on Bio-Rad iQ5 instrument). Melting analysis can be done immediately after the amplification reactions.

Number of Cycles	Step	Temperature	Time	Detection
1	Pre-denature	95 °C	10min	No
40	Denature	95 °C	10sec	No
	Annealing	See reference below	20sec	No
	Extension	72 °C	10sec	Yes

All-in-One™ miRNA qPCR Primer Validation Report


Temperature range	Rate of temperature change	Duration	Detection
66 °C ~ 95 °C	0.5 °C/ step	6 sec/ step	Yes
30 °C		30 sec	No

Note: The above conditions are designed for use with the Bio-Rad iQ5Q-PCR instrument. If other instruments are used, adjust the extension time and conditions for analysis as recommended by the manufacturers. The annealing conditions for each miRNA assay may be different. Please refer to the details of the conditions in the following section.


III. Test Results

1. hsa-miR-16 (hsmq-0031 primer)

Recommended annealing temperature : 60°C

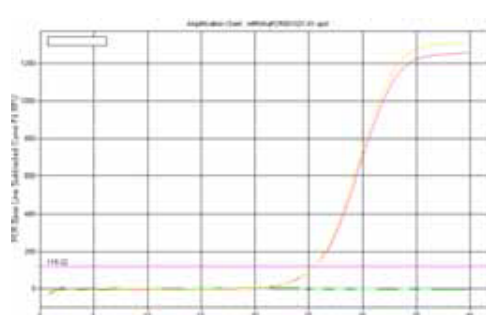


hsa-miR-16 Amplification Plot

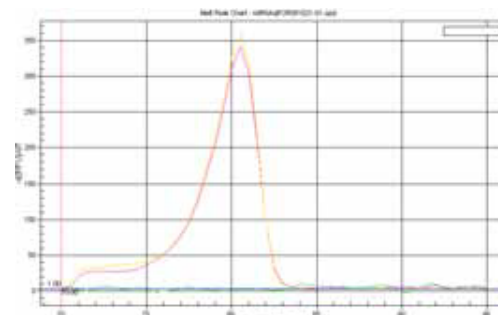


hsa-miR-16 melting analysis curve

2. hsa-miR-124 (hsmq-0032 primer)

Recommended annealing temperature : 60°C

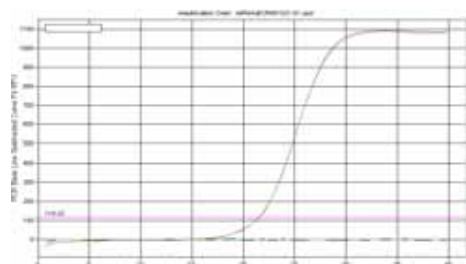

hsa-miR-124 Amplification Plot

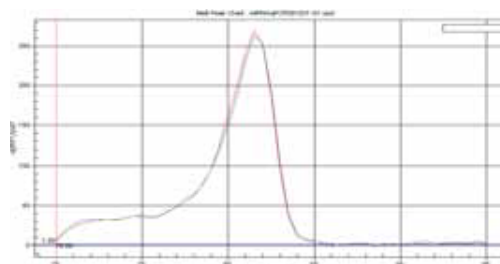


hsa-miR-124 melting analysis curve


All-in-One™ miRNA qPCR Primer Validation Report

3. hsa-miR-125b (hsmq-0034 primer)

Recommended annealing temperature : 60°C

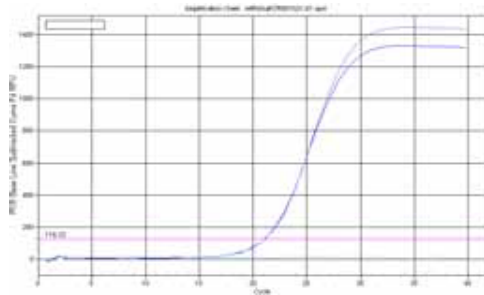


hsa-miR-125b Amplification Plot

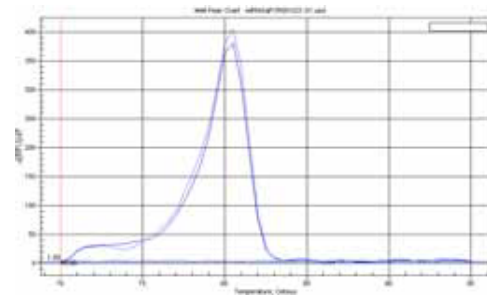


hsa-miR-125b melting analysis curve

4. hsa-miR-137 (hsmq-0035 primer)

Recommended annealing temperature : 60°C

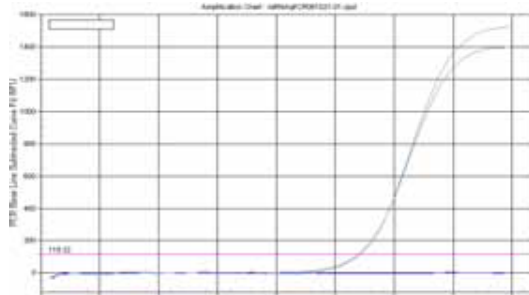


hsa-miR-137 Amplification Plot

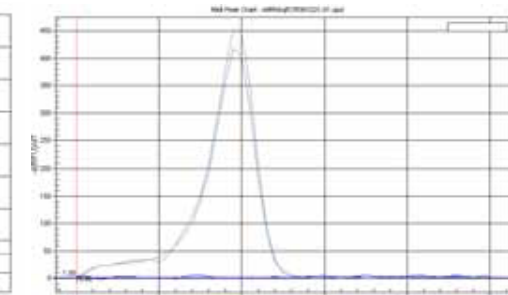


hsa-miR-137 melting analysis curve

5. hsnRNA U 6 primer testing

Recommended annealing temperature : 60°C

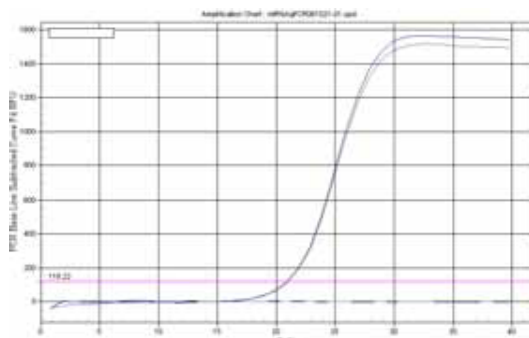


hsnRNA U6 Amplification Plot

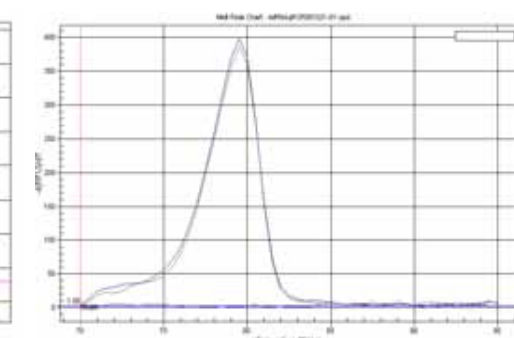

hsnRNA U6 melting analysis curve

All-in-One™ miRNA qPCR Primer Validation Report

6. Electrophoresis Results

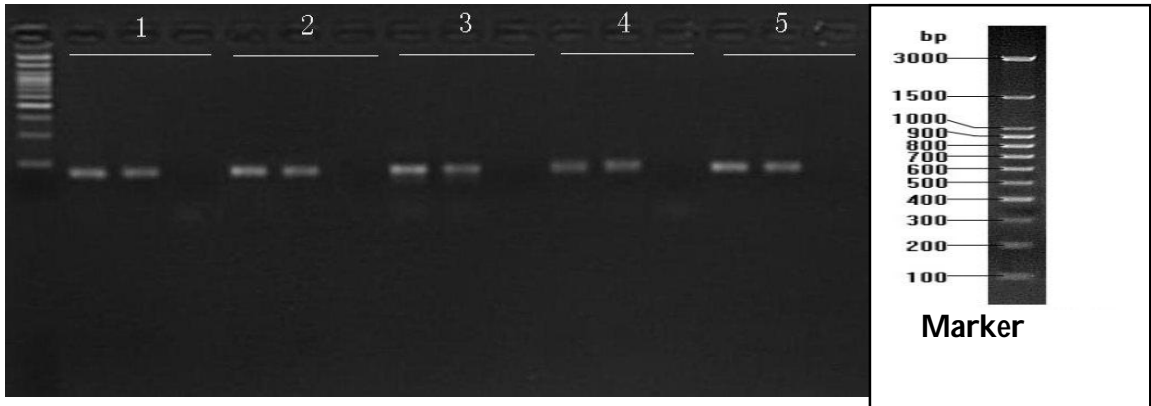


5 µl of the PCR of products, run on 3% agarose gel. Each set contains two positive controls and one NTC
 Lane 1: hsa-miR-16; Lane 2: hsa-miR-124; Lane 3: hsa-miR-125b; Lane 4: hsa-miR-137; Lane 5: hsnRNA-U6.

7. hsa-miR-19a (hsmq-0655 primer)


7.1 Validation.

The following link may be used to view a larger amplification plot and melting curve as well as a gel electrophoresis image of this primer's validation results. Contact GeneCopeia at inquiry@genecopeia.com or call 866-360-9531 to have an electronic document sent to you.

http://www.genecopeia.com/product/search/mirna_qpcr_validate.php?id=miRQP0293&mid=hsa-miR-19a

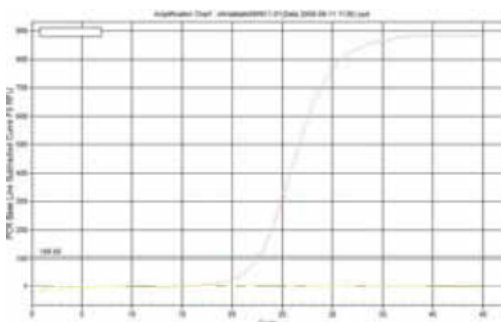

hsa-miR-16 Amplification Plot

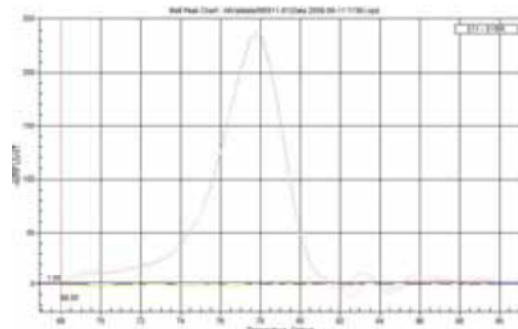

hsa-miR-16 melting analysis curve

7.2 Discrimination assay of two primers with single-base differences in their sequences


7.2.1 has-miR-19a BLAST

hsa-miR-19a	UGUGCAA <u>A</u> UCUAUGCAA <u>A</u> ACUGA
hsa-miR-19b	UGUGCAA <u>C</u> AUGCAA <u>A</u> ACUGA

The BLAST alignment shows that a single-base difference exists between the two miRNAs: hsa-miR-19a and miR-19b. To discriminate and validate the primer of interest (has-miR-19a) two plasmids were constructed for each of the two miRNAs.

All-in-One™ miRNA qPCR Primer Validation Report

7.2.2 Sequence result of the miRNA qPCR plasmids

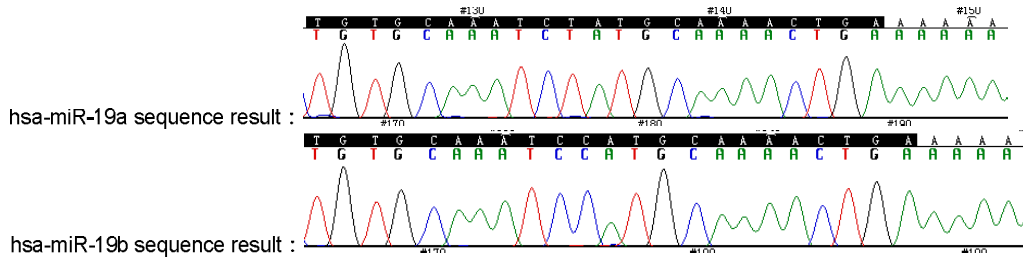

7.2.3 qPCR validation

Using the All-in-One miRNA qRT-PCR Detection Kit, the hsa-miR-19a-specific primer was validated using miRNA expression plasmids for has-miR-19a and has-miR-19b as templates (about 10^7 molecules/reaction).

The percent relative discrimination was calculated using the differences between the Ct values of the mismatched template and the matching template with respect to the hsa-miR-19a-specific primer (Relative discrimination = $2^{-\Delta Ct} \times 100\%$). The results are shown in the table below:

Annealing temperature °C	Ct value for miRNA expression plasmid amplification		Relative discrimination
	hsa-miR-19a	hsa-miR-19b	hsa-miR-19b
63.3	20.30	30.89	0.0651%
61.4	18.40	25.81	0.5880%
58.9	17.50	21.57	5.9545%

7.2.4 Conclusion

Based on these results, a qPCR annealing temperature between 59.9 and 61.4 °C is recommended. This range gives the highest amplification efficiency as well as the best relative discrimination between the two single-base-different miRNAs.

All-in-One™ miRNA qPCR Primer Validation Report

IV. Limited Use License and Warranty

Limited Use License

Following terms and conditions apply to use of all OmicsLink™ ORF Expression Clones in all lentiviral vectors and Packaging Kit (the Product). If the terms and conditions are not acceptable, the Product in its entirety must be returned to GeneCopoeia within 5 calendar days. A limited End-User license is granted to the purchaser of the Product. The Product shall be used by the purchaser for internal research purposes only. The Product is expressly not designed, intended, or warranted for use in humans or for therapeutic or diagnostic use. The Product must not be resold, repackaged or modified for resale, or used to manufacture commercial products without prior written consent from GeneCopoeia. This Product should be used in accordance with the NIH guidelines developed for recombinant DNA and genetic research. Use of any part of the Product constitutes acceptance of the above terms.

Limited Warranty

GeneCopoeia warrants that the Product meets the specifications described in the accompanying Product Datasheet. If it is proven to the satisfaction of GeneCopoeia that the Product fails to meet these specifications, GeneCopoeia will replace the Product. In the event a replacement cannot be provided, GeneCopoeia will provide the purchaser with a refund. This limited warranty shall not extend to anyone other than the original purchaser of the Product. Notice of nonconforming products must be made to GeneCopoeia within 30 days of receipt of the Product. GeneCopoeia's liability is expressly limited to replacement of Product or a refund limited to the actual purchase price. GeneCopoeia's liability does not extend to any damages arising from use or improper use of the Product, or losses associated with the use of additional materials or reagents. This limited warranty is the sole and exclusive warranty. GeneCopoeia does not provide any other warranties of any kind, expressed or implied, including the merchantability or fitness of the Product for a particular purpose.

GeneCopoeia is committed to providing our customers with high-quality products. If you should have any questions or concerns about any GeneCopoeia products, please contact us at 301-762-0888.

© 2009, GeneCopoeia, Inc.

GeneCopoeia, Inc.
9620 Medical Center Drive, #101
Rockville, Maryland 20850
Tel: 301-762-0888 Fax: 301-762-8333
Email: inquiry@genecopoeia.com
Web: www.genecopoeia.com

To order contact:

**306, Aggarwal City Mall, Opposite M2K Pitampura,
Delhi – 110034 (INDIA). Ph: +91-11-42208000, 42208111, 42208222
Mobile: +91-9810521400 Fax: +91-11-42208444
Email: customerservice@atzlabs.com
Web: www.atzlabs.com**