

RNA is stable in LifeGuard Solution

RNA Isolated from soil stored at -20° C, 4° C, and RT for 1 day, 1 week and 30 days in LifeGuard™

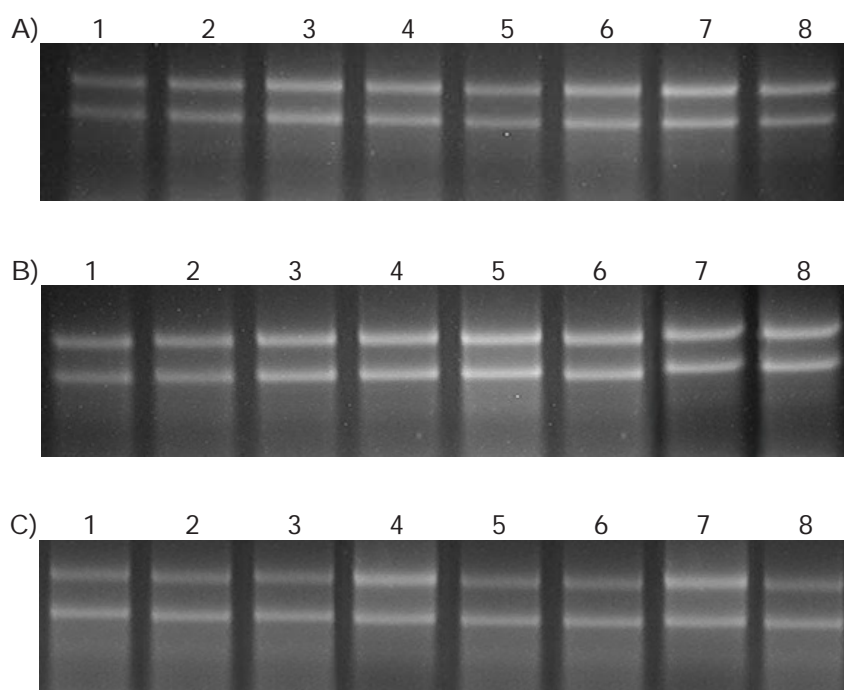


Figure 1 A-C A) Soils stored for 1 day in LifeGuard™ B) Soil stored in LifeGuard™ for 1 week, and C) soil stored for 30 days in LifeGuard™. All samples were extracted with the RNA PowerSoil® Kit. Numbers across the top indicate the storage temperature as follows: 1+2= -20° C, 3+4 = 4° C, 5+6 = room temperature, 7+8 = fresh soil stored at 4° C (control). detected.